OenoFoss Test Name Change


This procedure will walk you through the steps needed to change the name of any desired test.

Begin by navigating to Foss Integrator. Click the "Care" button, and then "Go to Foss Integrator"



Enter the password of "1" when prompted


Using the Menu bar, click "Window", and then "Settings

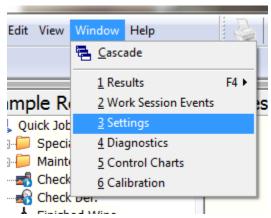


Double Click "Product Settings"




Expand the profile that contains the test you want to rename, and click "Prediction Models"

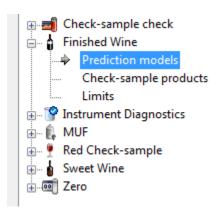

*Double click the name of the test you want to rename

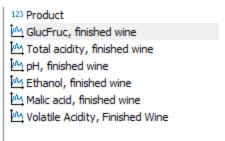

*Select the "Data Output" tab. The name listed under the "Header" column is the display name for the test. Change it to your desired name.

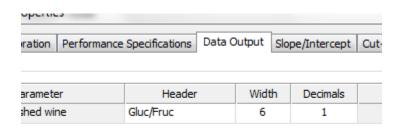

orce negatives to zero

*Press OK when done.

Repeat the steps marked with a (*) to rename any additional tests.

Press OK on the Product Settings windows when all changes are complete.

OK

End